

PRODUCT INFORMATION

Decalcifier concentrate plus indicator

Cat. No.: 18141

PRODUCT DESCRIPTION

Decalcifier is a pH neutral reagent to remove calcium from tissue specimen on the basis of a complex former.

Usage:

Working dilution: 1:2 (1+1) to 1:10 (1+9)

The working dilution depends on the calcium amount that should be dissolved. The Decalcifier has a calcium-specific indicator. If it turns from red to blue, it will be necessary to add more Decalcifier solution.

Storage: at +15 °C to +30 °C

Vers. 09/09