

Protease-Inhibitor-Mix M

Cat.No. : 39102

Cat. number	Components	Target Protease
12745	AEBSF-HCl	Serine Proteases
13718	Aprotinin from bovine lung	Serine Proteases
14980	Bestatin	Aminopeptidase B and Leucine Aminopeptidase
21100	E-64	Cysteine Proteases
51867	Leupeptin	Cysteine Proteases and Trypsin-like Serine Proteases
52682	Pepstatin A	Aspartic Proteases
Storage (°C)		-15 to -25
Reconstitution	Dissolve the content of the vial (15.8mg) in 1ml DMSO to obtain a 100x stock solution (suitable for the treatment of 100ml tissue extract). After reconstitution, aliquot and freeze at -20°C (stable for at least 4 weeks).	

Die Eignung des Produktes für spezielle Anwendungszwecke wird nicht zugesichert.
We do not guarantee that the product can be used for a special application.