

---

**Protease-Inhibitor-Mix FY**

**Cat.No. : 39104**

---

<b>Cat. number</b>	<b>Components</b>	<b>Target Protease</b>
12745	AEBSF-HCl	Serine Proteases
21100	E-64	Cysteine Proteases
52682	Pepstatin A	Aspartic Proteases
32010	1,10-Phenanthroline	Metalloproteases
<b>Storage (°C)</b>		-15 to -25
<b>Reconstitution</b>	Dissolve the content of the vial (62.6mg) in 1ml DMSO to obtain a 100x stock solution (suitable for the treatment of 100ml tissue extract). After reconstitution, aliquot and freeze at -20°C (stable for at least 4 weeks).	

**Die Eignung des Produktes für spezielle Anwendungszwecke wird nicht zugesichert.**  
We do not guarantee that the product can be used for a special application.