

Protein Test Mixture 6

Cat.No. : 39207

Protein Components	Molecular weight
Phosphorylase B	97 400
Albumin bovine (BSA)	67 000
Albumin egg	45 000
Carbonic anhydrase	29 000
Trypsin inhibitor (soybean)	21 000
Cytochrome C	12 500
Trypsin inhibitor (bovine lung)	6 500
Storage (°C)	+2 to +8

Reconstitution Dissolve to a concentration of 1mg/ml in SDS-sample buffer.
A volume of 5µl is sufficient for SDS-PAGE (in a standard form of 80 x 60 x 0.75mm).

SDS-sample buffer:

125 mM Tris-HCl pH 6.8; 2% SDS; 15% glycerol;
10 mM DTT; 0.025% Bromphenol blue; 0.025% Orange G

Die Eignung des Produktes für spezielle Anwendungszwecke wird nicht zugesichert.
We do not guarantee that the product can be used for a special application.