

Protein Test Mixture 5 for SDS PAGE

Cat.No. : 39209

Protein Components	Molecular weight (kDa)
Trypsin inhibitor bovine lung	6 500
Cytochrom C	12 500
Trypsin inhibitor soybean	21 000
Carbonic anhydrase	29 000
Storage (°C)	+2 to +8

Reconstitution

Dissolve to a concentration of 1mg/ml in SDS-sample buffer.
A volume of 5µl is sufficient for SDS-PAGE (in a standard
form of 80 x 60 x 0.75mm).

SDS-sample buffer:

125 mM Tris-HCl pH 6.8; 2% SDS; 15% glycerol;
10 mM DTT; 0.025% Bromphenol blue; 0.025% Orange G

Die Eignung des Produktes für spezielle Anwendungszwecke wird nicht zugesichert.
We do not guarantee that the product can be used for a special application.