

Product of the Month


SERVA
■ serving scientists ■

Get **20%** discount* now!

April

*For all SERVA customers within the EU. Outside Germany and Austria online shop orders only. Can not be combined with other discounts, promotions etc.

SingleQuant Assay Kit Fast and Reliable Protein Quantification

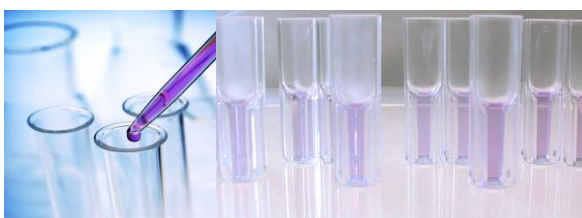

SingleQuant Assay Kit provides precise and reproducible analysis data of protein quantification in single tube format. Reducing agents, e.g. DTT, 2-Mercaptoethanol or detergents, e.g. SDS, CHAPS do not interfere with the assay method.

Features:

- Suitable for **2 – 1400 µg protein**
- Completed in only 45 min
- Stable absorption for min. 2 hours

The assay method bases on the pH-dependent precipitation of proteins with acidic, ethanolic amido black 10B solution (Popov et al., *Acta Biol. Med. Ger.* 1975: 34(9): 1441 – 1446).

Ordering information

Product	Cat. no.	Size	List price	Special price*
SingleQuant Assay Kit	39226.01	200 Tests	159.00 €	127.20 €

* The special price is valid until April, 30th 2021.

For further information, please contact:

Dr. Judith Koch
Tel.: +49 6221 13840-44
E-Mail: Tech.Service@serva.de

Workflow:

5 min	Precipitation of protein-dye complexes
	↓
5 min	Centrifugation
	↓
10 min	Washing of protein-dye complexes, 2x
	↓
15 min	Resolubilization of the dye with NaOH
	↓
10 min	Measurement of absorption at 624 nm

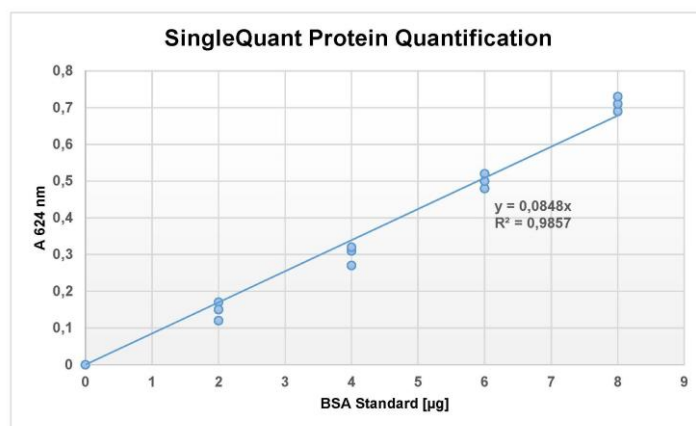


Fig.: Calibration curve with bovine serum albumin (BSA) standards

SERVA Electrophoresis GmbH
Carl-Benz-Str. 7
D-69115 Heidelberg • Germany
www.serva.de